



Inline Particle Sizing



Inline Particle Sizing for Process Control of Fluid Bed and High Shear Mixing Processes

Dr. Michael Koehler, Parsum GmbH Chemnitz

17th International Drying Symposium 2010

Wigratec Session

October 3 – 6, 2010, Magdeburg





Presentation Outline


- Terms of On-Line, In-line, At-line etc.
- Destination and Requirements
- Presentation of Particle sizing results
- PARSUM's measurement principle
- Case Study Fluid Bed Processes
- Case Study High shear granulation processes
- Conclusion





In-line Particle Size Analysis (PSD) Definition of Terms

- **On-line** – Measurement with “real time” output of the results (e.g. by continuous or quasi- continuous sampling or bypass)
- **In-line** – Measurement and analysis performed directly inside the process
- **AT-line** – Measurement of a sample near the process in production environment
- **Off-line** – All laboratory methods with sampling (analysis is separated from process in respect to time and space)

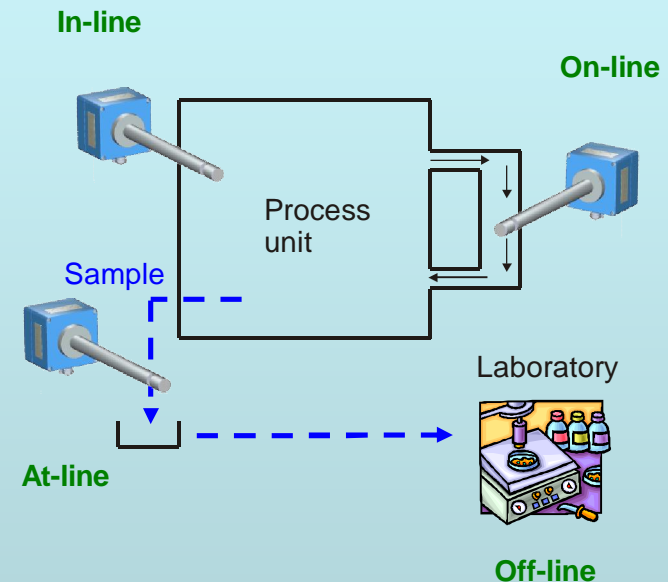



Why In-line Particle Size Analysis ?

- Monitors the particle size continuously
- Gives more representative results
- Increases efficiency and throughput.
- Gives higher quality by tighter control
- Reduces the potential product contact to operator
- Improves batch to batch consistency
- Gives better control to minimize wastage
- Reduces costs of production





Requirements for use in Industry


- Low hardware efforts for installation
- Wide measuring range ... μm – ... mm
- True inline capability
- Robust industrial design
- Long term stability
- Independent of process conditions and product properties



Presentation of Particle Sizing Results

Off-line (process photo):

- PSD only at sample times
- No information about the progress of particle size

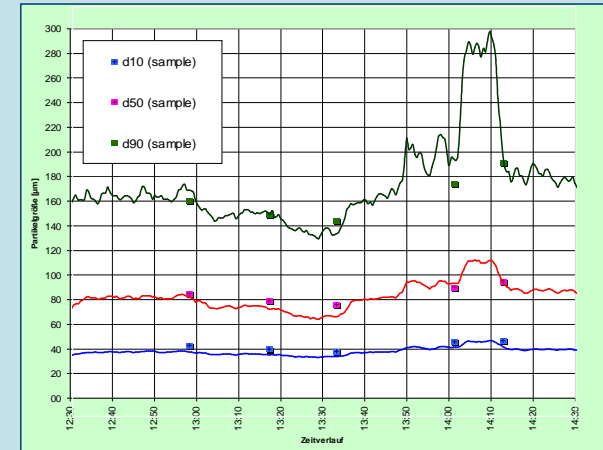



In-line (process movie):

- Progress of the PSD over the time
- Concentrated information for process control e.g. $X_{10,3}$; $X_{50,3}$ and $X_{90,3}$
- Dynamic of measurement is adjustable to process dynamic
- Averaged PSD at any interval





Parsum measuring principle

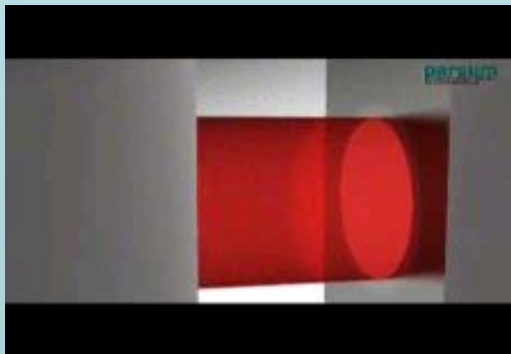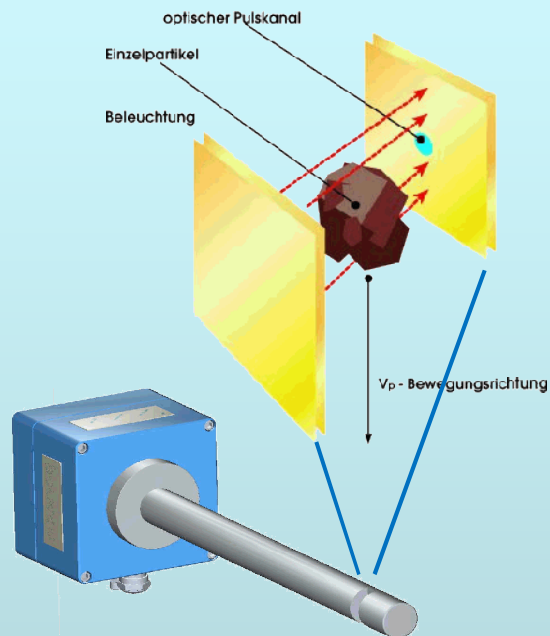


- Modified spatial filter method
- Measuring principle is based on evaluation of the shadows from moved particles in a parallel laser beam

Properties:

- Measuring range approx. 50µm to 6mm
- Velocity range up to 50 m/s
- Continuously recording of particle velocity and particle size
- Method especially for online use
- Chord length measurement
- Single particle measurement
- Probe with no moving parts



Modified Spatial Filter Method

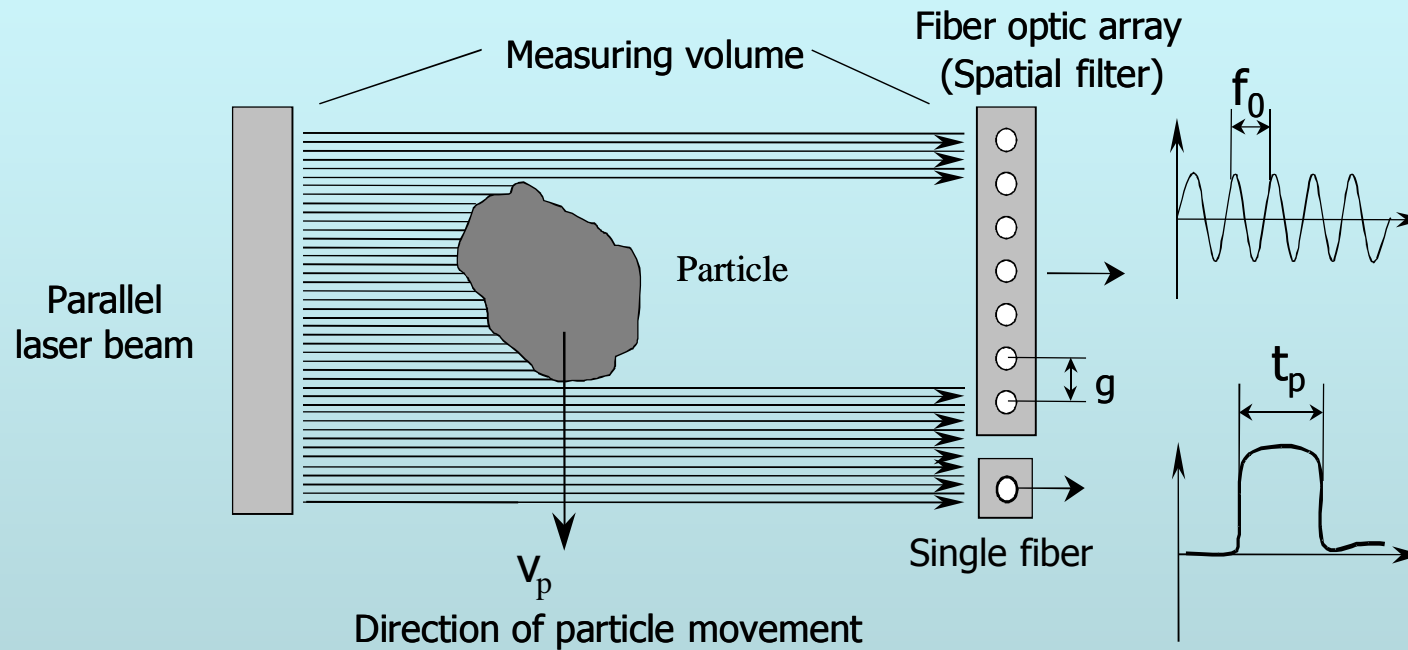

$$\text{Velocity: } v_p = f_0 * g$$

$$\text{Size: } x_p = t_p * v_p - d$$



Process Interface – Inline Eductor D23



- For high loaded particle streams and for processes with irregular particle movement like Fluid Bed, High Shear Granulation
- Air purge to hold optics free
- Ring-nozzles to suck and accelerate particles
- Particles are moving on straight Lines
- separation effect is minimized
- Installation with Tri Clamp DN50
- Periodic back purge function with a separate internal nozzle



Fluid Bed Batch Granulation


Objective:

- To see trends in granulation,
 - To find end-point of granulation,
 - To see any defect (blocked nozzles, break down of fluidized bed ...)
and
 - To demo the usability of IPP-70 for direct measurement inside fluidized beds.
-
- Product: Lactose Powder,
 - Equipment: 5 Kg lab scale FB Granulator, Top Spray
 - Installation: IPP70-S with D23






Fluid Bed Batch Granulation

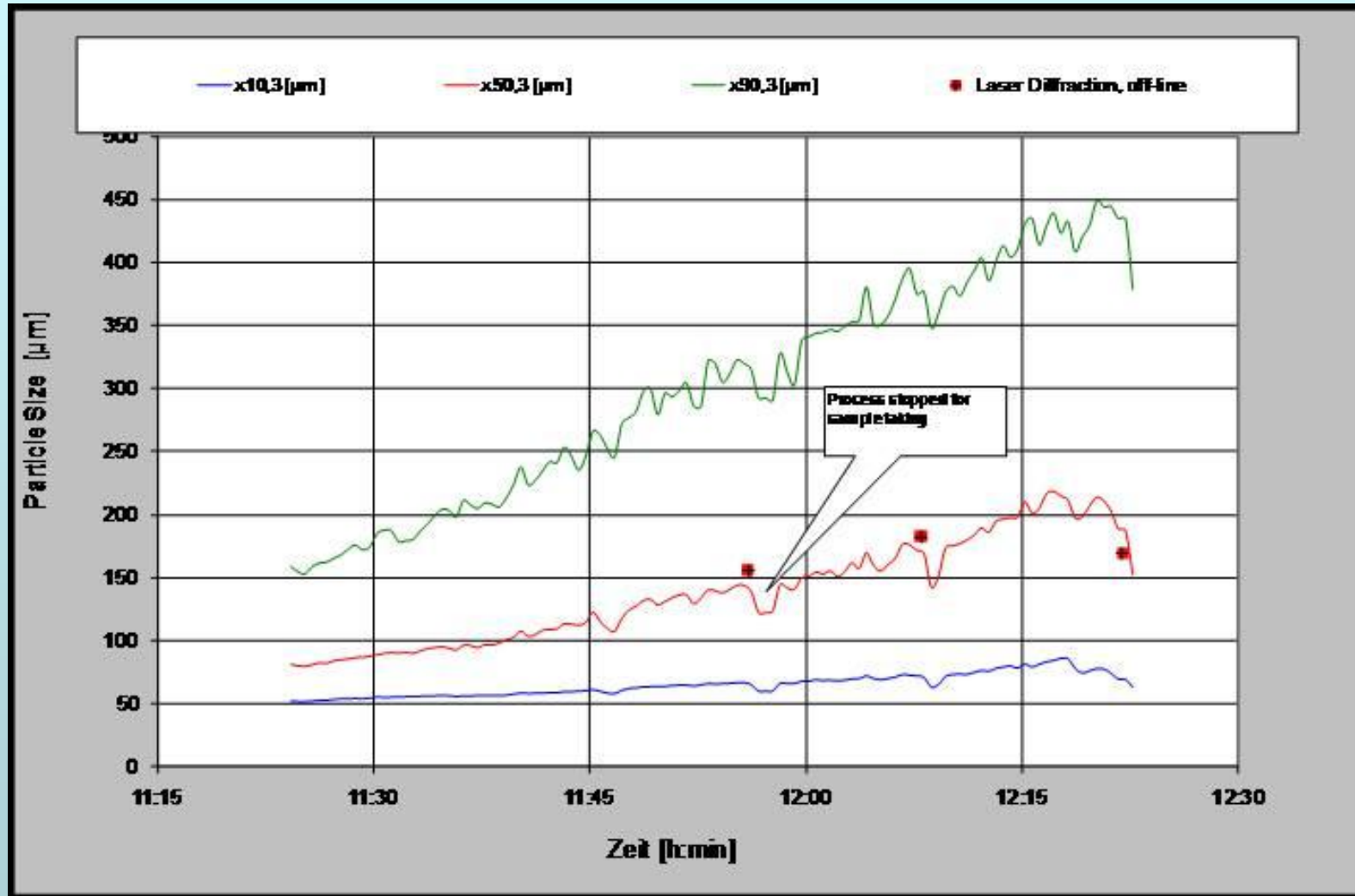



Wurster Coating

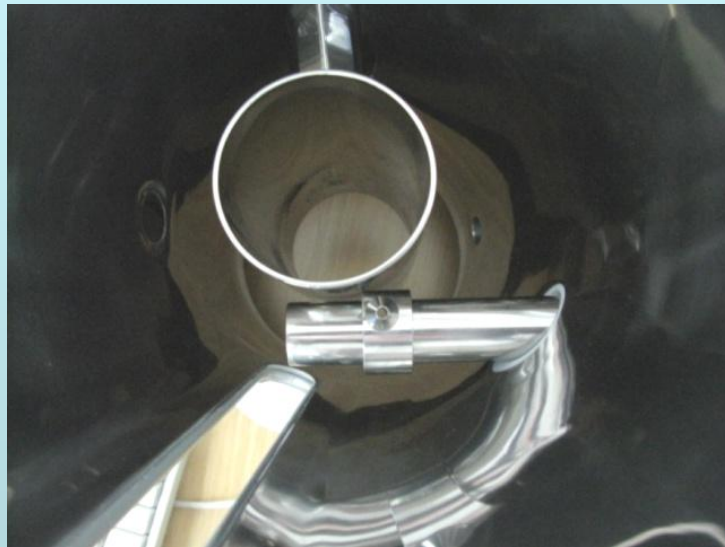

Objective: Measurement of thickness of sprayed layer
Detection of Agglomerates (Twins)


Product: Sugar pellets, spherical, 300µm,

Process-Equipment: Lab-scale FB-Granulator, 3Kg with Wurster-Tube and Bottom-Spray

Installation: IPP70-S with Inline-eductor D23




Wurster Coating

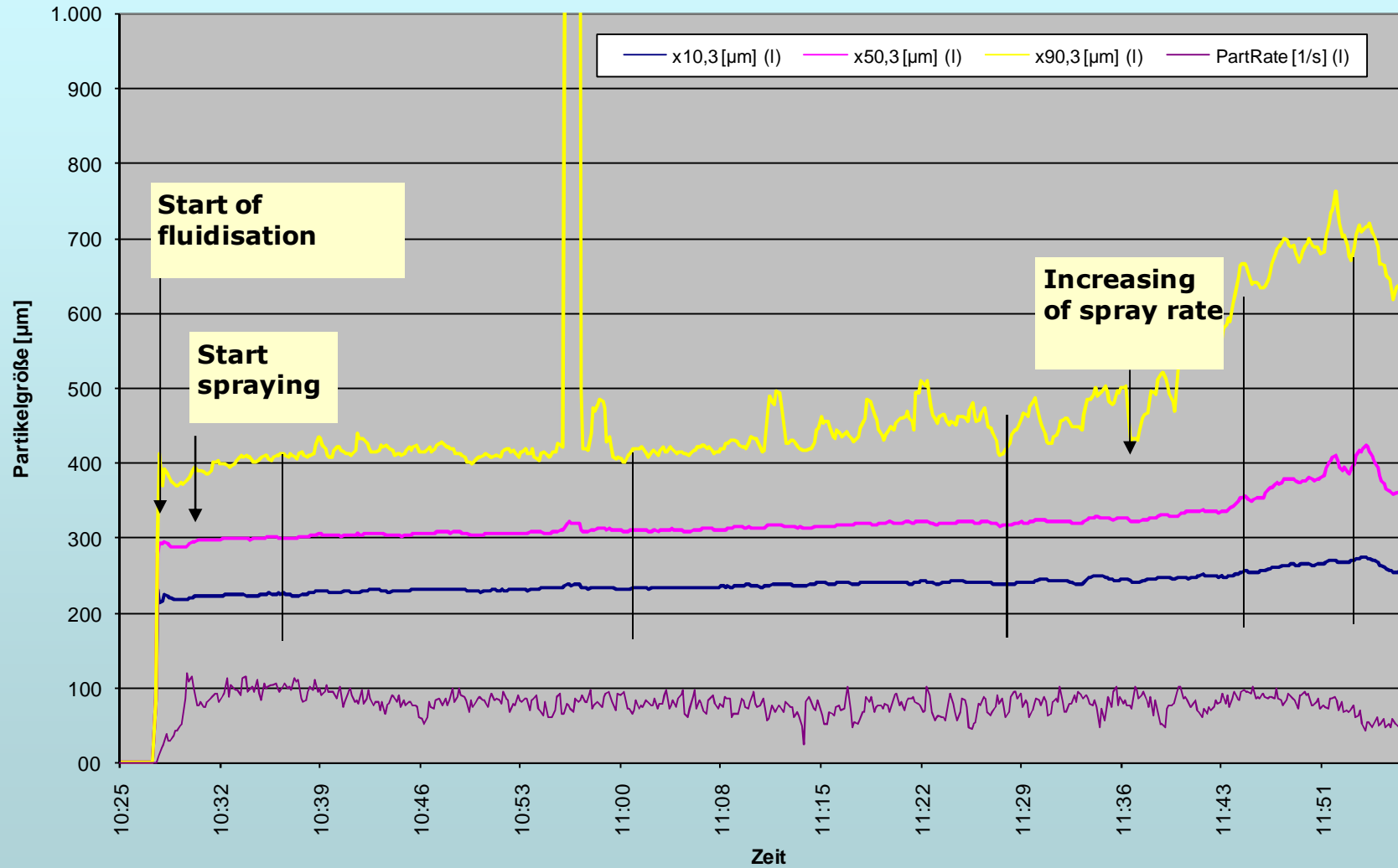



Wurster Coating

— q3 density dist. 10:36:36 — q3 density dist. 11:01:16 — q3 density dist. 11:28:36 — q3 density dist. 11:45:16 — q3 density dist. 11:53:06

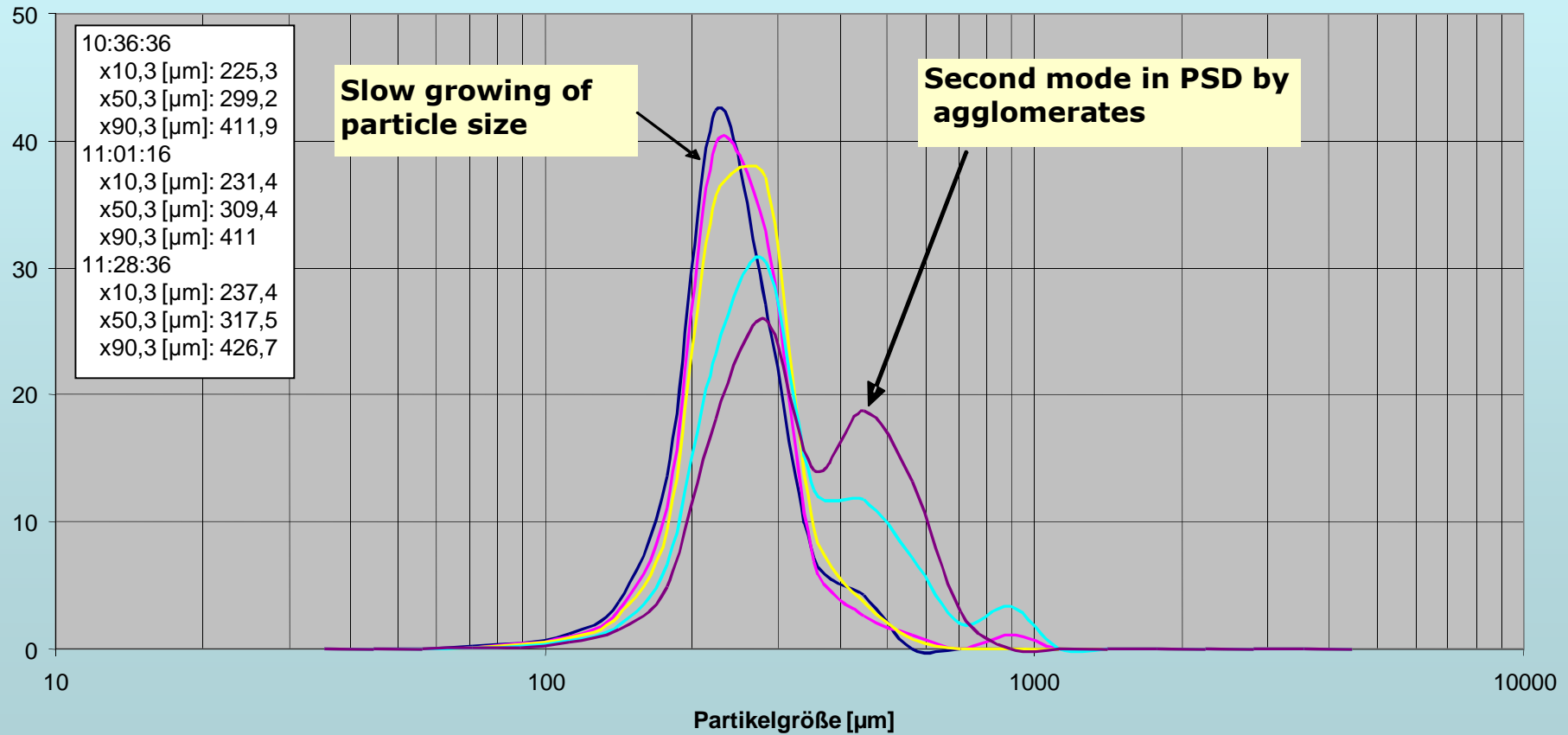



High Shear Granulation, production scale



Objective:

To demo useability of IPP
70-SL in larger mixers.
To follow fast particle size
changing
flexible depth of probe

Product:

Lactose with API


Installation:

IPP 70-SL (60 cm)
with inline-eductor D23






High Shear Granulation, production scale

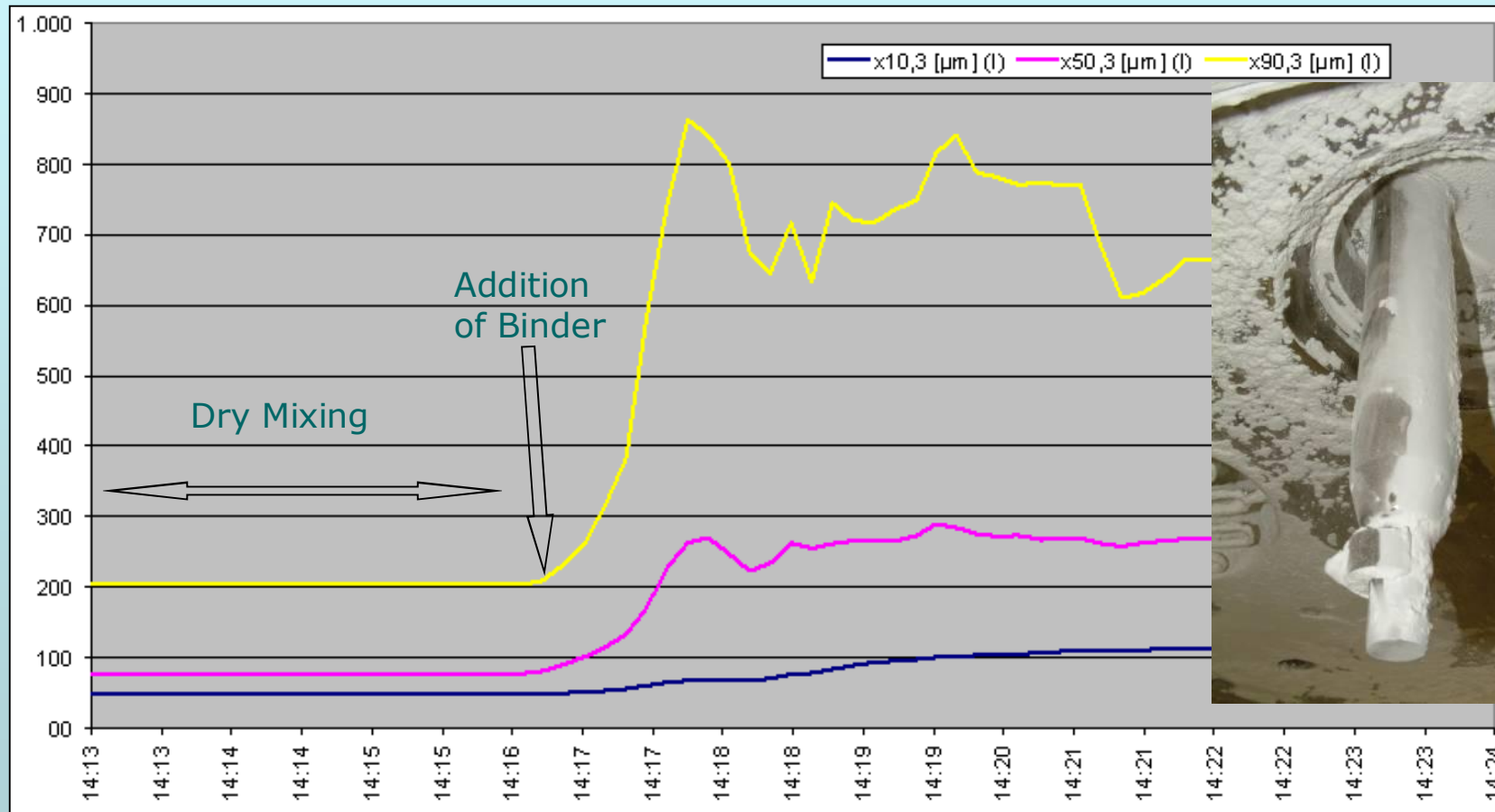



High Shear Granulation, production scale

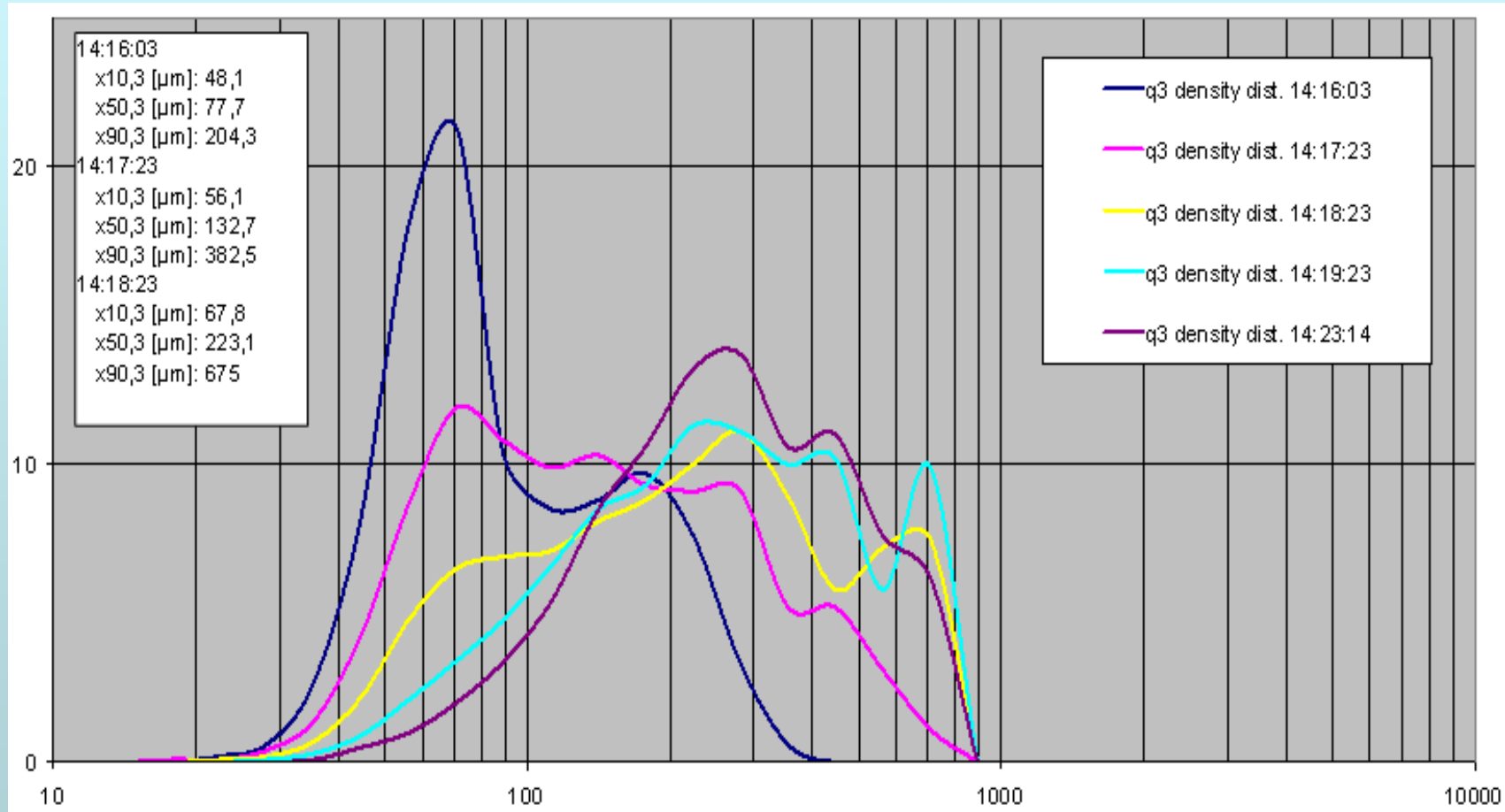



References

- D. Petrak, Simultaneous Measurement of Particle Size and Particle Velocity by the Spatial Filtering Technique, *Part. Part. Syst. Charact.*, 19, 391-400 (2002).
- S. Schmidt-Lehr, H.-U. Moritz and K. C. Jürgens, Online Control of Particle Size during Fluidised Bed Granulation, *Pharm. Ind.*, 69, 478-484 (2007).
- T. Närvänen, T. Lipsanen, O. Antikainen, H. Räikkönen and J. Yliruusi, Controlling granule size by granulation liquid feed pulsing, *Int J Pharm.*, 357, 132-138 (2008).
- A. Burggraeve, T. Van Den Kerkhof, M. Hellings, J. P. Remon, C. Vervaet and T. De Beer, Evaluation of in-line spatial filter velocimetry as pat monitoring tool for particle growth during fluid bed granulation, *Eur. J. Pharm. Biopharm.*, doi: 10.1016/j.ejpb. 2010.06.001.
- J. Huang, C. Goolcharran, J. Utz, P. Hernandez-Abad, K. Ghosh and A. Nagi, A PAT Approach to Enhance Process Understanding of Fluid Bed Granulation Using In-line Particle Size Characterization and Multivariate Analysis, *J Pharm Innov*, 5, 58-68 (2010) DOI 10.1007/s12247-010-9079-x.
- Jarvis Jeejing Ng, In-line particle size analysis for high shear granulation processes, AFACT 10, 28.04.2010 – 30.04.2010, Manchester
- D. Petrak, S. Dietrich, G. Eckardt and M. Köhler, In-line particle sizing for process control by an optical probe based on the spatial filtering technique (SFT), in: CD Proc. 6th World Congress Particle Technol., H H 1 0 00613, Nuremberg (2010).